


Fetal Bovine Serum

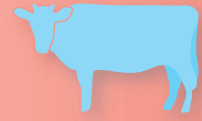

- High Quality, Reliable FBS.**
- Free Lot Check Samples.**
- Free Storage.**
We offer a free storage for your FBS.
We can ship them as you need.
- Customer can use same batch of FBS in US and Japan.**



Nichirei has continued distributing FBS for over 30 years in Japan. Our FBS will powerfully support your international research!

USDA approved FBS *1

	Product name	Product Code	volume	Price
NICHIREI	FBS	174012-500ML	500mL	quote

*1 Please contact us about country of origin and price. We offer several countries of origin for FBS.

Nichirei Biosciences is a member of International Serum Industry Association.



INTERNATIONAL SERUM INDUSTRY ASSOCIATION



Product Properties

Filtration	Triple 0.1 μ m sterile filter
Storage conditions	-10 to -40 ° C
pH	6.8 – 8.1
Osmolality	260 -330 mOsm/kg H ₂ O
Sterility (Bacteria & Fungi)	Not Detected
Mycoplasma	Not Detected
Hemoglobin	< 25 mg/dL
Endotoxin	< 100 EU/mL
Bovine IgG	Record
Tetracycline	Record
Total Protein	3.0 -4.5 g/dL
Protein gel electrophoresis	Record
Growth Promotion	Tested
Virus(9CFR)	Record
Biochemical prof.	Record



Safety FBS

- ◆ Raw materials are collected from healthy animals.
- ◆ Raw materials are tested and confirmed to meet product criteria.
- ◆ Products are filtered through sterile filter 3 times.



Support

- ◆ **Recommended Thawing method**
 1. Leave the bottles at room temperature for 10 minutes.
 2. Set the water bath to 30-37C in advance and put the bottles in it.
 3. Swing the bottles slightly every 10-15 minutes and thaw the FBS completely.
- ◆ **Heat inactivated (HI) FBS**
 Recommended HI condition is 56C \pm 1C for 30 minutes.
 (We can provide HI FBS.
 Please note that Lot check sample is non-HI.)

Please feel free to contact us.

NICHIREI BIOSCIENCES INC.

6-19-20, Tsukiji, Chuo-ku Tokyo, 104-8402, Japan
 Phone: +81-3-3248-2207
 Fax: +81-3-3248-2243
 E-mail: n1060x006@nichirei.co.jp
 Site: <https://www.nichirei.co.jp/bio/english/>

[Distributor]

Nacalai USA, Inc.

10225 Barnes Canyon Rd., Suite A103 San Diego, CA 92121, USA
 Phone: (858) 404-0403
 Fax: (858) 404-0408
 E-mail: info@nacalaiusa.com
 Site: <https://www.nacalaiusa.com/>