
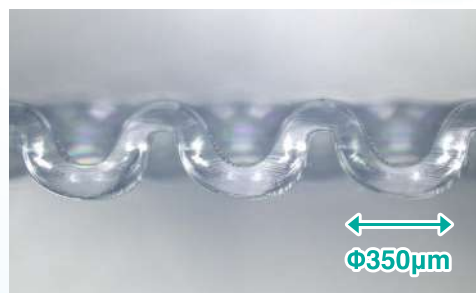


WELLBAG-S350

Now
on
sale

WELLBAG-S350 has 32,000 microwells with a well diameter of 350 μ m. The size of the bag is the same with WELLBAG-S500, so the same dedicated holder can be used. Since this product has the same well diameter with a large size WELLBAG: WELLBAG-L350, it can be used for scaleup studies.

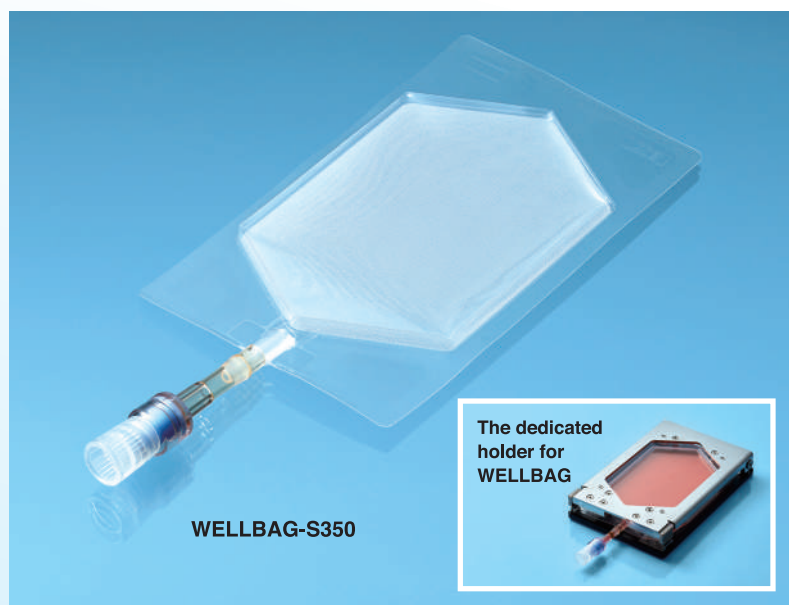

Features

- ① Low contamination risk due to the closed system
- ② Mass culture of uniform size spheroids
- ③ Efficient recovery of formed spheroids
- ④ Easy gas exchange during culture by using highly gas permeable film
- ⑤ Spheroid migration can be prevented by using a dedicated holder
 - ※The size of the bag is the same with WELLBAG-S500, so the same dedicated holder can be used.
- ⑥ Safety standards according to the Japanese Pharmacopoeia
- ⑦ To remove air bubbles in the wells, incubation method is recommended.
 - ※Please refer the operation manual of WELLBAG for details.

For the Research and Development,
Manufacturing of Cell Therapeutics

Closed
system

A large number
of uniform size
Spheroids

Easy recovery
of spheroids



Toyo Seikan Group Holdings, Ltd.

Comparison of WELLBAG-S350 and WELLBAG-S500

	WELLBAG-S350	WELLBAG-S500
Size	72 mm×120 mm	
Material	Polyethylene with non-cell adhesion coating	
Culture Area	50 cm ²	50 cm ²
Culture volume	10~20 mL	10~20 mL
Well diameter	350 μm	500 μm
Number of wells	ca.32,000	ca.18,000




※ WELLBAG-S350 is stamped with 32K on the seal of the bag.

Suitable for the scale up study of spheroid culture

- The large size WELLBAG: WELLBAG-L350 has a culture area of 1,000cm² and 650,000 wells with a diameter of 350μm. Standard medium volume is 400ml.
- WELLBAG-S350 has a culture area of 50cm² and 32,000 wells with a diameter of 350μm. The standard medium volume is 20ml, which is approximately the same amount of medium per well as the large size WELLBAG.

Air bubble removal for WELLBAG-S350

- Incubation method is recommended. It takes about a half day to remove air bubbles by this method.
 - By placing this product filled with the specified amount of culture medium in a special holder and leaving it overnight in an incubator set at 37°C. Air bubbles in the wells will naturally be released to the outside of the bag.
 - Seed the cell suspension into the bag from which air bubbles have been removed.
- ※ Because the well diameter of S350 is smaller than that of S500, it is difficult to remove air bubbles manually.

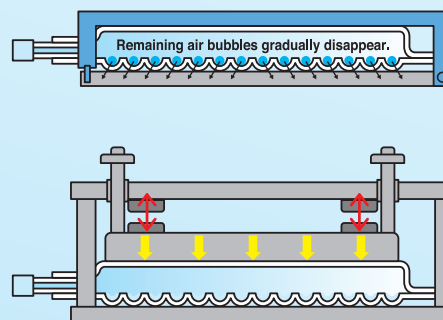

Note : The use of WELLBAG to insulin-producing cells including pancreatic precursor cells may be restricted depending on the intended use. So, please contact the manufacturer or a distributor in advance for this use.



Toyo Seikan Group Holdings, Ltd.

Business Management Group, Business Innovation Office
2-18-1 Higashi-Gotanda, Shinagawa-ku, Tokyo, Japan
Tel : +81-3-4514-2305 Fax : +81-3-3280-8121
(MAIL) bis_innov_wellbag@tskg-hd.com
(URL) <https://www.tskg-hd.com/cellsolut/>

※ The contents are subject to change without notice. Please note this in advance.

Dealer :

Nacalai USA, Inc.

6625 Top Gun Street, Suite 107, San Diego, CA 92121
TEL: (858) 404 0403
Email: info@nacalaiusa.com
Website: www.nacalaiusa.com